
SST, SST Science

Tables

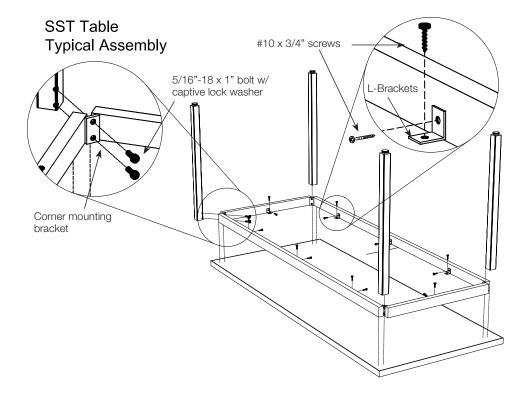

Hardware Included:



L-Bracket



#10 x 3/4" screws (in to rails and wooden tops)



#8 x 3/8" screws (used to attach frame to epoxy tops)



5/16"-18 x 1" PH bolt w/captive lock washer

Tools Required:

- #2 Phillips Head Screwdriver
- #3 Phillips Head Screwdriver

Step 1:

Attach each leg to the apron by bolting through a corner mounting bracket into welded threaded insert in leg using (2) 5/16-18 x 1" bolt per leg.

Step 2:

To attach frame assembly to the table top place the table top upside down on a clean, padded surface. Center the length and width of welded table frame on the underside of the top. Secure frame to table top through L Brackets with #10 x 3/4" PH Screws into both the frame and bottom of top.

